

~~TOC~~

## Diamond NV Centers

### Basics

see also [W Nitrogen-vacancy center](#)

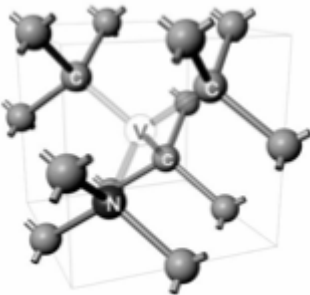


Figure1: Model of the nitrogene vacancy in the diamond lattice

### Excitation

- wavelength 532nm
- excitation power should be 100  $\mu$ W or above

### Emission

Figure 3 shows the emission spectra of a NV-center. The pronounced Peak at 637nm is the zero-phonon line of the diamond lattice (comparable to Raman).

- Emission filters therefor should be 650LP.

### Fluorescence Lifetime

- Fluorescence Lifetime is approx. 12ns.

## Measuring NV centers

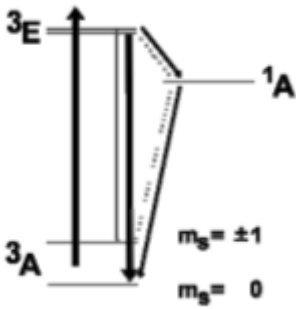


Figure2: Jablonski diagram of the luminescent transition of the nitrogen-vacancy in a diamond lattice. Bright fluorescence corresponds to the  ${}^3E \rightarrow {}^3A$  ( $m_s=0$ ) transition

Due to their photo-physical stability the NV-centers make for excellent demonstration samples for [antibunching measurements](#). They are also potential candidates to image the [confocal volume](#).

The following will explain how to find and measure NV-centers in bulk diamond.

## System setup

- refractive index of diamond is 2.4. that means oil immersion is necessary as well as a high N.A.
- used objective: N.A. 1.3 is working, better 1.45 oil!
- Although measurements were possible also with the water immersion objective: with  $60\times 1.2$  water less than 50% count rate in comparison to  $100\times 1.3$  oil
- 0.95 N.A. air objective did not yield any usable results.
- excitation source: 532 nm
  - MOFA (40MHz - for finding the centers, later 10MHz (12ns fluorescence lifetime))
- excitation power: >7000 a.u. @ 40MHz, > 2000 a.u. @ 10MHz (basically as high as possible)
- major dichroic: 532 / 635
- emission bandpass: 690/70 (for [antibunching measurements](#) 2 bandpasses directly in front of detectors)
- objective: 1.3 N.A. oil immersion

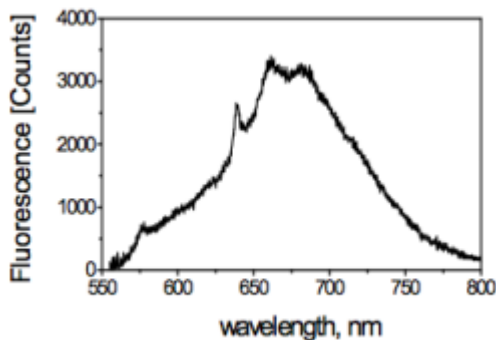


Figure3: fluorescence spectrum of a single NV center in diamond excited with 532 nm

## Mounting the sample

sample is mounted directly on top of the objective (with immersion oil, without any coverslips etc.).

## Imaging/ finding nv centers

- focus on the interface between oil and diamond (nice reflection should be visible due to refractive index mismatch between oil and diamond)
  - when imaging the surface most likely fluorescent spots will be found. Those are probably dirt and unlike the NV-centers will bleach in a matter of seconds. They also have a shorter fluorescence lifetime.

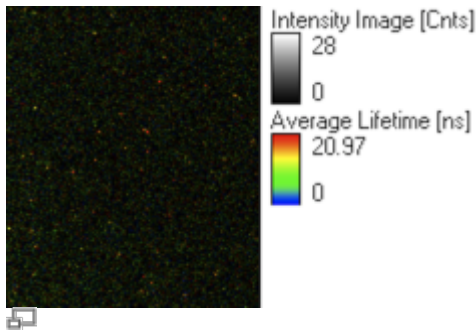


Figure4: Prescan of NV-centers in bulk diamond. 5 micrometer into the sample, 80×80 micrometer, 150×150 pixel resolution. 532nm excitation, 690/70 emission, 40MHz rep. rate, approx. 7000 a.u.

- move approx. 5 micrometer into the sample.
- Although photo-physical stability makes the nv-centers ideal single emitter samples compared to single molecules the laser power to excite them has to be considerably higher. (0.1 mW and above).
- Figure 4 shows a prescan of the complete field of view. The orange dots are NV-centers.
- zoom into a region with a few NV-centers (see figure 5),
- select one, and position the focus on the NV-center.
- use the oscilloscope and pifoc+piezo to move the NV-center in x,y and z into the center of the focus.
- For lifetime analysis switch laser repetition rate to 10MHz (see figure 6)
- check whether the fluorescence lifetime is around 12 ns.
- the NV-center's emission should stay stable (no bleaching)

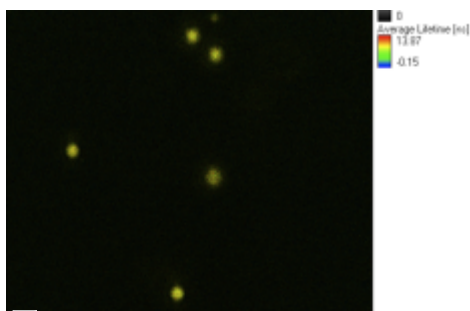


Figure5: NV-centers in bulk diamond. 5 micrometer into the sample, 532nm excitation, 690/70 emission, 40MHz rep. rate, 6789.96a.u., 2ms/pixel, 15×15 micron

## Typical Measurement Results

- Prescan should look like figure 4
- smaller area is depicted in figure 5 and figure 6
- Selecting a ROI in figure 6 and plotting the fluorescence lifetime histogram should resemble figure 7
- when measuring the fluorescence intensity of a single NV center, the count rate should stay stable. The time trace in figure 8 shows a reduction of the countrate due to stage drift (temperature etc, or in this case probably because the MT200 inhouse II was mounted on a wheelbarrow instead of an optical table)
- to perform [antibunching measurements](#) on a selected NV-center choose a NV-center that is isolated, measuring time of 120 s should be enough to yield an antibunching measurement comparable to figure 9.

*NOTE:* Measurement setup has to be changed (see [antibunching measurements](#))

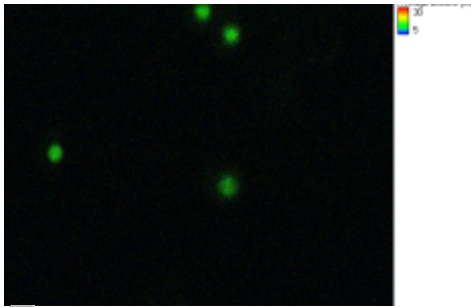


Figure6: same as figure 5 but recorded with 10 MHz rep rate., 2304.94 a.u.;4 ms/pixel 45 min rec. time!

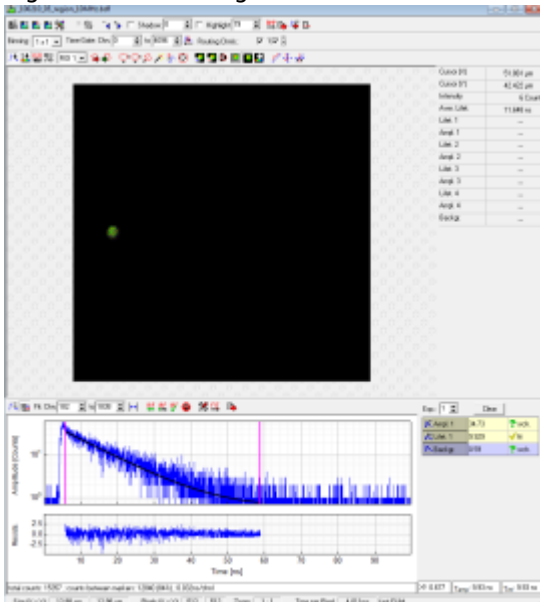


Figure7: Fluorescence Lifetime Fit of a selected NV center, 10 Mhz rep rate, Inhouse II, MPD/SPAD

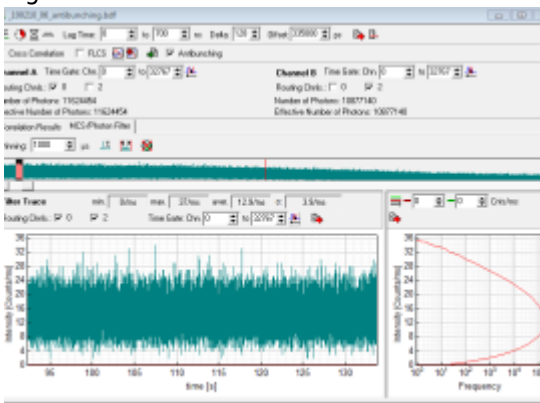


Figure8: No bleaching visible. reduction in count rate is due to mechanical drift (inhouse II) on wheelbarrow

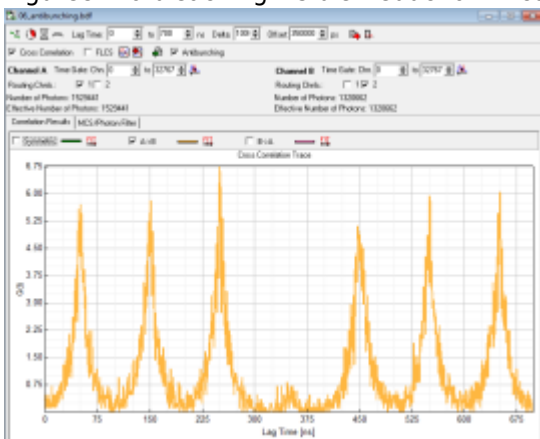


Figure9: Antibunching measured on NV-center. Acquisition time 120 s, 1.3 N.A. oil immersion, 3534.98 a.u. 532

## References

M. Boersch, R. Reutera, G. Balasubramaniana, R. Erdmann, F. Jelezkoa, J. Wrachtrup Fluorescent nanodiamonds for FRET-based monitoring of a single biological nanomotor FoF1-ATP synthase **Proc. of SPIE Vol. 7183, 71832N** (2009)

<http://dx.doi.org/10.1117/12.812720>

Copyright of this document belongs to PicoQuant GmbH. No parts of it may be reproduced, translated or transferred to third parties without written permission of PicoQuant GmbH. All information given here is reliable to our best knowledge. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications and external appearances are subject to change without notice.



PicoQuant GmbH  
Rudower Chaussee 29 (IGZ)  
12489 Berlin  
Germany

P +49-(0)30-1208820-89  
F +49-(0)30-1208820-90  
info@picoquant.com  
www.picoquant.com